

MWM EXTERNAL PROBE – HOW TO CHANGE THE BATTERY?



1. Remove **BOTH** end caps off the white tube as follows:







2. Remove / slide the circuit board from inside the white tube (DO NOT disconnect the sensor that is attached):



3. Take out the old battery (take note of which way the battery is to be inserted - Positive points to the wires). Now fit the NEW Battery.





- 4. Slide the circuit board back into the white tube and replace / tighten both end caps.
- 5. Place your MWM Sensor back into your Fridge / Freezer now
- 6. LOGIN to the MATOS Monitoring website https://matosmonitoring.com/ and check you can see your fridge / freezer temperatures coming through live!