



MATOS[®] WIRELESS MONITORING KIT (MWM)

MATOS MWM Wireless Tag Setup Instructions

*Your MWM kit has already been preconfigured and assigned to the MATOS Monitoring Webserver.
In order to start temperature monitoring, please follow the below steps.*

To start tag monitoring:

1. Remove tag manager from packaging and plug in USB Power Cable.
2. Plug in the tag manager to an Ethernet/Network connection.
(Note: the Tag Manager operates in DHCP mode, so no additional configuration is required).
3. Once the 'cloud' connection indicator symbol on the front of the tag manager has turned solid green, the Tag Manager has established a connection with the web server.
4. Pull the plastic tab from the MWM Wireless Tag to begin monitoring automatically.
5. Place the tag inside the refrigerator or asset to be monitored. Please ensure it is in range of the Tag Manager (within 100m or less is advised).

To view/download temperature information:

1. Open a web browser and navigate to matosmonitoring.com.
2. Login with your account credentials
3. You should now see live temperature information on the Overview or Dashboard pages.
4. To check/configure alert contact information, click on the "Alert Contacts" Link.
5. To view historical temperature data, click the "Reports" link.

***You should now see all your live temperature data on
the MATOS Monitoring Website!***



Need Assistance:

Please do not hesitate to contact your Sales representative and/or the Service Team via emailing service@rollexmedical.com or by calling 1300 880 441 if you have any questions or need assistance.